Pharmacy Refrigerator DIN 13277 Advance Line

Modelo · Model

MPRA 66 G

Ref.

800902V1400

EAN: 5601359902141

HSCode: 84185019

W 450
D 470
H 740

60

38

+2/+8

Descrição

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Title Pres.

Equipment developed for use in pharmacy, for the conservation of drugs and vaccines, in accordance with normative DIN 13277. Due its dimensions, it can be placed on top of a counter or suspended in the wall using the metal bars on the back. The thermoformed plastic interior facilitates the cleaning, preventing the accumulation of bacteria and germs, and the exterior made of special coated steel prevents the appearance of corrosion with use.

Excellent stability, uniformity and temperature accuracy, guaranteed through the 50 mm water-based polyurethane insulation, refrigeration system with ventilated cold and automatic defrosting, as well as a quick temperature recovery after opening the door. The factory setpoint is +5°C, which can easily be adjusted in the range between +2°C and +8°C. With internal temperature homogeneity of +/- 2°C. It has a hole with diameter 12mm to pass a probe for external monitoring system.

Standard with lock to prevent unauthorized access, 3 removable shelves in plastic coated steel to prevent corrosion and basket on bottom, reversible glass door with low emissive double tempered glass and no condensation system, with automatic return to the closed position, led lighting, handle with ergonomic design and easy cleaning, and 4 adjustable feet allowing to level the equipment.

Características Advance line

Termostato Advance Advance Controller

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with keyboard lock.

Audible and visible alarms

High temperature
Low temperature
Open door
Probe failure
Power failure

Alarm Test

Test of the functionality of the alarms

Safety controller

Double safety to prevent the temperature from dropping from +2°C

Real time clock

Year, month, hour, minutes and seconds for data recording definition.

Equipment identification

Definition of the equipment serial number in the thermostat for register control.

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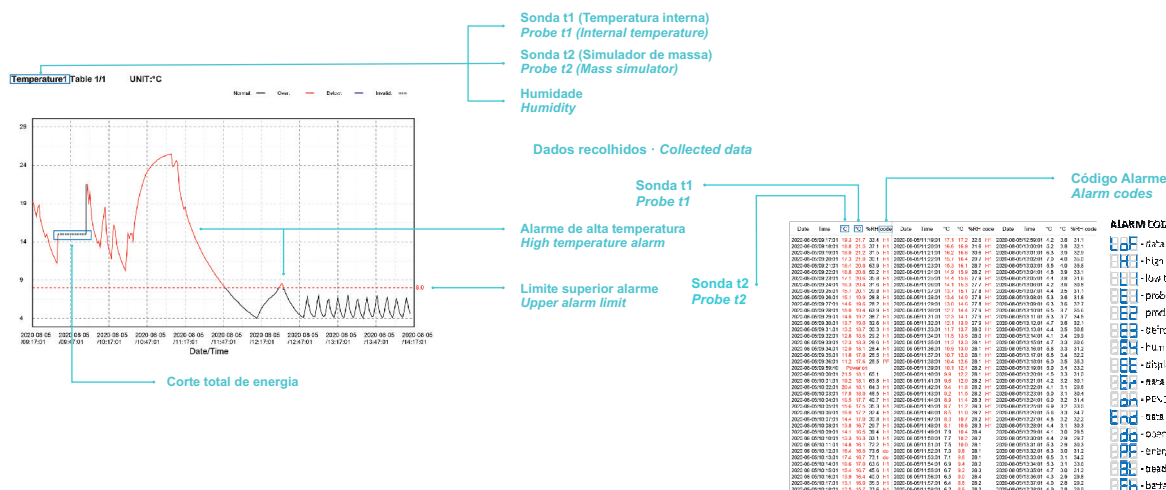
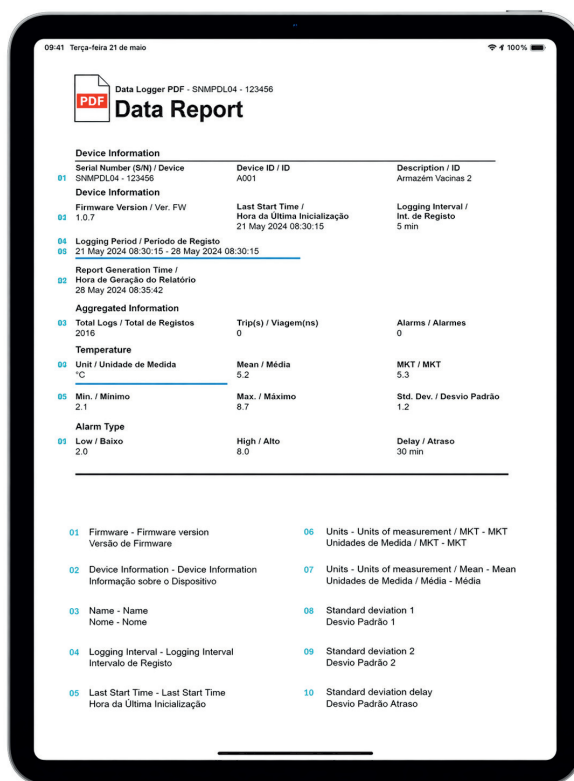
The Advance thermostat records the unit's temperature performance at two measurement points, as well as alarm events, with readings taken every 10 minutes and storage capacity for up to two years of data. These records can be downloaded via the USB port, generating a PDF file as shown in the example below.

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