Pharmacy Refrigerator DIN 13277 Advance Line

Modelo · Model

MPRA 350 S

Ref.

800903V1400

EAN: 5601359903148

HSCode: 84185019

W 600
D 700
H 1510

354

86

+2/+8

Descrição

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medgree.com

Title Pres.

Equipment developed for use in pharmacy, for the conservation of drugs and vaccines, in accordance with normative DIN 13277.

Excellent stability, uniformity and temperature accuracy, guaranteed through the 50 mm water-based polyurethane insulation, refrigeration system with ventilated cold and automatic defrosting, as well as a quick temperature recovery after opening the door. The factory setpoint is +5°C, which can easily be adjusted in the range between +2°C and +8°C. With internal temperature homogeneity of +/- 2°C. It has a hole with diameter 12mm topass aprobe for external monitoring system. The interior and exterior in special plasticized steel prevents the appearance of corrosion with use.

Standard with lock to prevent unauthorized access, 5 removable shelves in plastic coated steel to prevent corrosion and basket on bottom, reversible solid door with automatic return to the closed position, led lighting, handle with ergonomic design and easy cleaning, and 4 rollers with 2 levellers for easy equipment movement.

A proud member of  Olitrem family

Características

Advance line

Termostato Advance Advance Controller

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with keyboard lock.

Audible and visible alarms

High temperature
Low temperature
Open door
Probe failure
Power failure

Alarm Test

Test of the functionality of the alarms

Safety controller

Double safety to prevent the temperature from dropping from +2°C

Real time clock

Year, month, hour, minutes and seconds for data recording definition.

Equipment identification

Definition of the equipment serial number in the thermostat for register control.

Humidity control (OPTIONAL)

If active, it allows monitoring the humidity in the inside the equipment

Datalogger

Query maximum/minium temperatures and last alarms in display.

Product probe

Probe with mass simulator, allowing to simulate the product temperature in any part of the equipment.

USB

Records are downloaded as PDF file, thus ensuring that the data cannot be changed.

Records

Graphical and tabular record of the following data: Probe 1 Temperature; Temperature Probe 2 (mass simulator); Alarms; Humidity (if you activate the function). *with readings every 10 minutes and with a capacity for 2 years of data (adjustable until 10 years with readings each 60 minutes).

Battery

In case of power failure. Capable of keeping alarms and logs active for up to 26 hours.

Remote alarm output and RS 485

Tension free contact for connection to external alarms and RS 485 for connection to building management systems

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