# Pharmacy and Laboratory Refrigerator

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[medgree.com](http://medgree.com)

## Title Pres.

Versatile equipment that can be used for both pharmacy and laboratory.

Excellent stability, uniformity and temperature accuracy, guaranteed through the 80 mm water-based polyurethane insulation, refrigeration system with air forced ventilated cold and automatic defrosting, as well as a quick temperature recovery after opening the door. The factory setpoint is +5°C, which can easily be adjusted in the range between +2°C and +16°C. With internal temperature homogeneity of +/-2.5°C. It has a hole with diameter 15mm to pass a probe for external monitoring system. The interior in stainless steel and exterior in special plasticized steel prevents the appearance of corrosion with use.

Standard with lock to prevent unauthorized access, 8 removable shelves (4 each door) in plastic coated steel to prevent corrosion and separators in the middle between shelves to the product don't fall, double tempered low-emission glass door with automatic return to the closed position, led lighting, handle with ergonomic design and easy cleaning, and 6 swivel (2 with brake) for easy equipment movement.

A proud member of  Olitrem family

# Características

## Advance line

### Termostato Advance Advance Controller

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with keyboard lock.

##### Audible and visible alarms

High temperature  
Low temperature  
Open door  
Probe failure  
Power failure

##### Alarm Test

Test of the functionality of the alarms

##### Safety controller

Double safety to prevent the temperature from dropping from +2°C

##### Real time clock

Year, month, hour, minutes and seconds for data recording definition.

##### Equipment identification

Definition of the equipment serial number in the thermostat for register control.

##### Humidity control (OPTIONAL)

If active, it allows monitoring the humidity in the inside the equipment

##### Datalogger

Query maximum/minium temperatures and last alarms in display.

##### Product probe

Probe with mass simulator, allowing to simulate the product temperature in any part of the equipment.

##### USB

Records are downloaded as PDF file, thus ensuring that the data cannot be changed.

##### Records

Graphical and tabular record of the following data: Probe 1 Temperature; Temperature Probe 2 (mass simulator); Alarms; Humidity (if you activate the function). \*with readings every 10 minutes and with a capacity for 2 years of data (adjustable until 10 years with readings each 60 minutes)

##### Battery

In case of power failure. Capable of keeping alarms and logs active for up to 26 hours.

##### Remote alarm output and RS 485

Tension free contact for connection to external alarms and RS 485 for connection to building management systems

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

The Advance thermostat records the unit's temperature performance at two measurement points, as well as alarm events, with readings taken every 10 minutes and storage capacity for up to two years of data. These records can be downloaded via the USB port, generating a PDF file as shown in the example below.

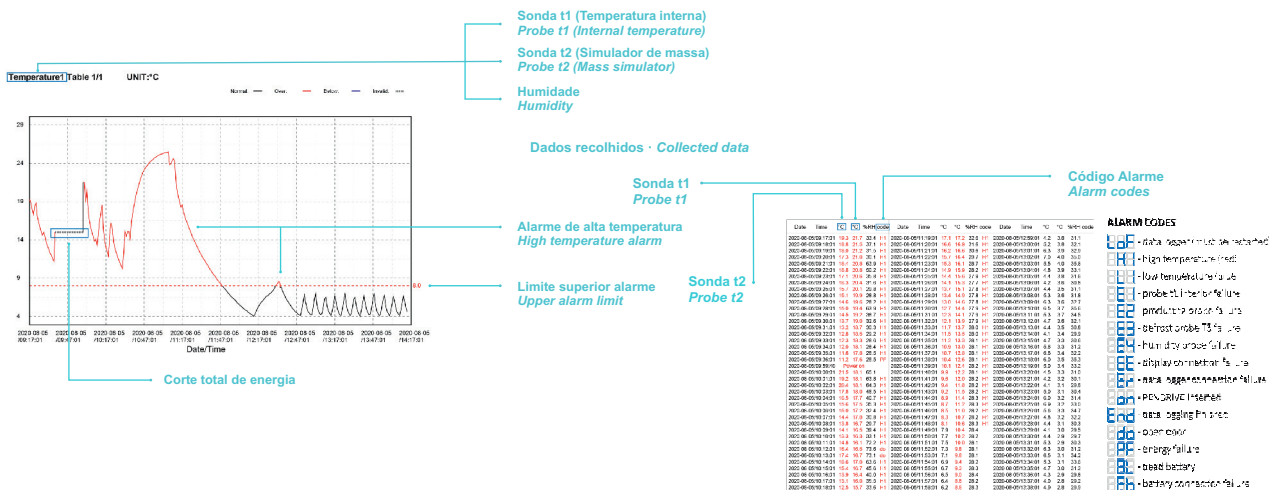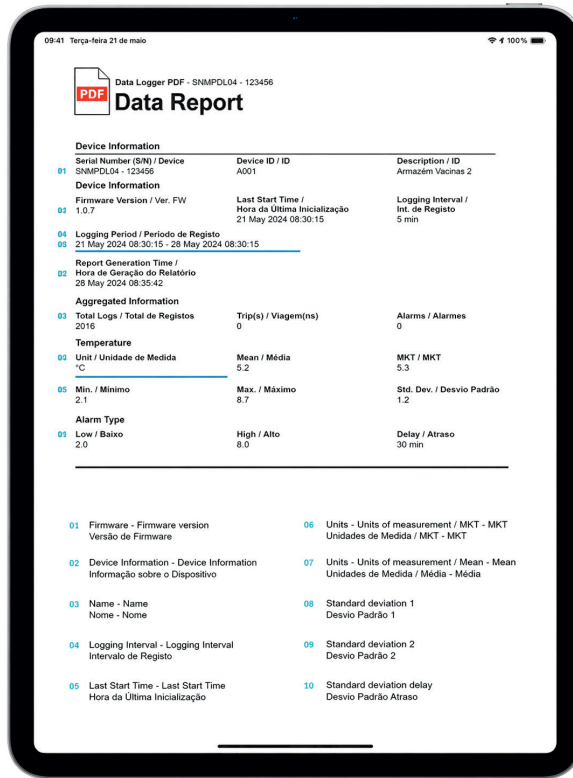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# Pharmacy and Laboratory Refrigerator Advance Line


Modelo · Model

MLRA 1400 G

Ref.

801083V100


## Accessórios · Accessories

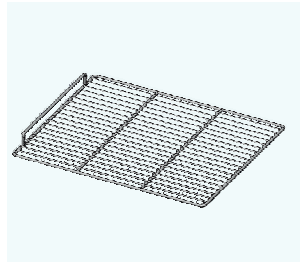
Prateleira Shelf  
Cód. 521455\_008


Kit Calha Suporte Prateleira · Shelf Holder Kit  
Cód. 129038\_63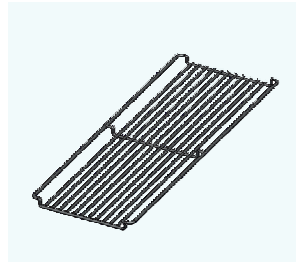

Separador de Prateleira · Shelf Separator  
Cód. 018009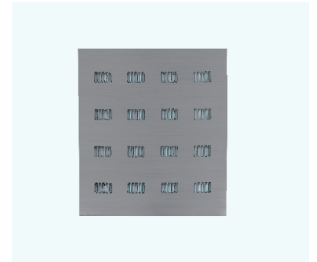

Prateleira Inox · Inox-Shelf  
Cód. 107464

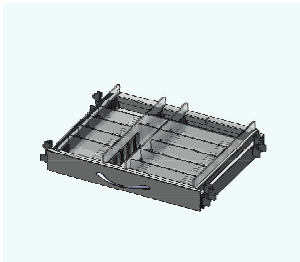Kit Gaveta Drawers Kit  
Cód. 129043V0749

Exterior em Inox Outside Inox  
Cód. 129045V0729

Embalagem Madeira Wood Package

# Pharmacy and Laboratory Refrigerator Advance Line

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