Laboratory Freezer Advance Line

Modelo · Model

MLFA 700 S

Ref.

801087V100

EAN: 5601359087107

HSCode: 84185011

W 720
D 840
H 2040

535

133

-10/-25

Descrição

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medgree.com

Title Pres.

Equipment developed to store samples in perfect safety conditions in laboratories, hospital and clinics.

Excellent stability, uniformity and temperature accuracy, guaranteed through the 80 mm water-based polyurethane insulation, refrigeration system with air forced ventilated cold and automatic defrosting, as well as a quick temperature recovery after opening the door. The factory setpoint is -18°C, which can easily be adjusted in the range between -10°C and -25°C. With internal temperature homogeneity of +/-5°C. The interior in stainless steel AISI 304 and exterior in plasticized steel prevents the appearance of corrosion with use. It has a hole with diameter 15mm topass aprobe for external monitoring system.

Standard with lock to prevent unauthorized access, 4 removable shelves in plastic coated steel to prevent corrosion, reversible solid door with automatic return to the closed position, led lighting when opening door, handle with ergonomic design and easy cleaning, and 4 swivel (2 with brake) for easy equipment movement.

A proud member of  Olitrem family

Características Advance line

Termostato Advance Advance Controller

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with keyboard lock.

Audible and visible alarms

High temperature
Low temperature
Open door
Probe failure
Power failure

Alarm Test

Test of the functionality of the alarms

Real time clock

Year, month, hour, minutes and seconds for data recording definition

Equipment identification

Definition of the equipment serial number in the thermostat for register control

Humidity control (OPTIONAL)

If active, it allows monitoring the humidity in the inside the equipment

Datalogger

Query maximum/minium temperatures and last alarms in display

Product probe

Probe with mass simulator, allowing to simulate the product temperature in any part of the equipment

USB

Records are downloaded as PDF file, thus ensuring that the data cannot be changed

Records

Graphical and tabular record of the following data: Probe 1 Temperature; Temperature Probe 2 (mass simulator); Alarms; Humidity (if you activate the function). *with readings every 10 minutes and with a capacity for 2 years of data (adjustable until 10 years with readings each 60 minutes)

Battery

In case of power failure. Capable of keeping alarms and logs active for up to 26 hours

Remote alarm output and RS 485

Tension free contact for connection to external alarms and RS 485 for connection to building management systems

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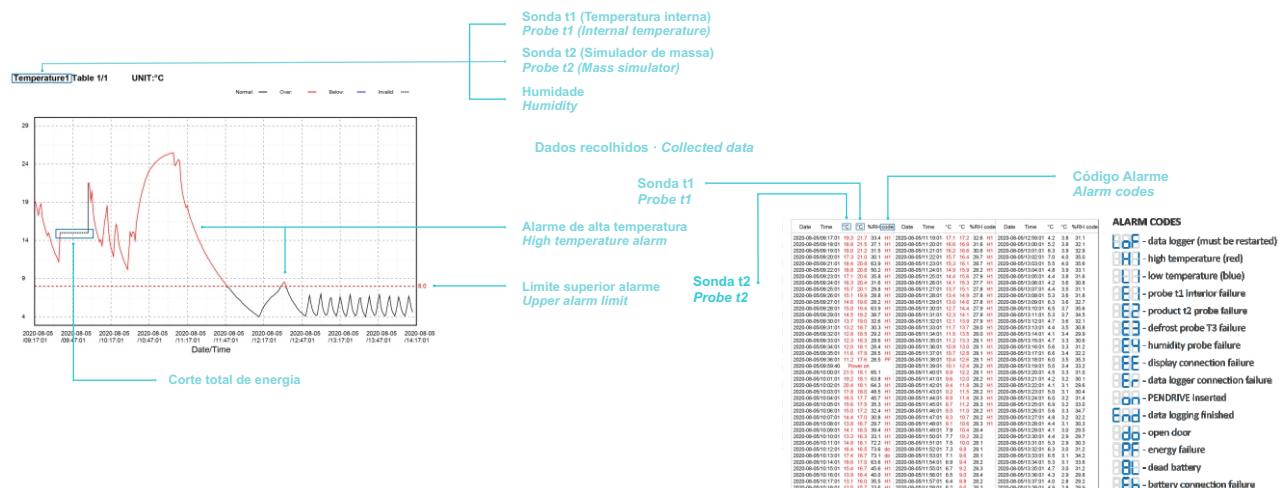
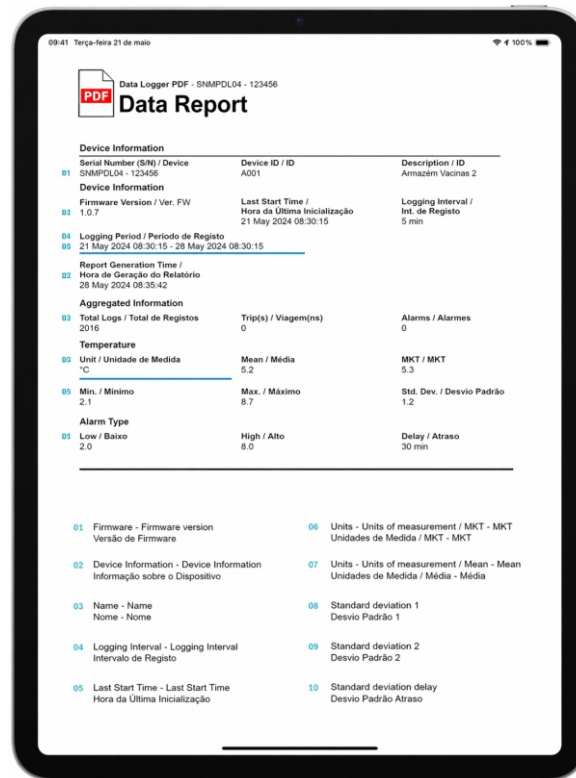
The Advance thermostat records the unit's temperature performance at two measurement points, as well as alarm events, with readings taken every 10 minutes and storage capacity for up to two years of data. These records can be downloaded via the USB port, generating a PDF file as shown in the example below.

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